

## Begin WiFi Setup

6




On device Press/  
HOLD WiFi button for  
3 secs.  
“AP” will appear  
on display.

Wait for a few seconds  
then press (no need to  
hold) WiFi button,  
“AP READY”  
will appear on display.


## Connect Unit to WiFi network via Mobile device

7

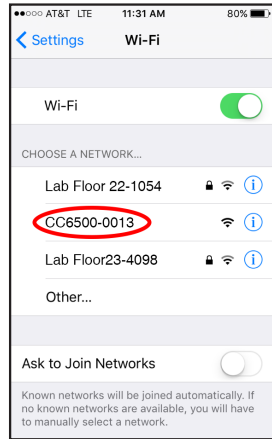


Using a WiFi enabled  
device (smart phone,  
tablet, etc.), find  
**CC65XX-XXXX**  
in available networks.

XXXX will be  
last 4 digits of unit's  
S/N# (on back of unit).

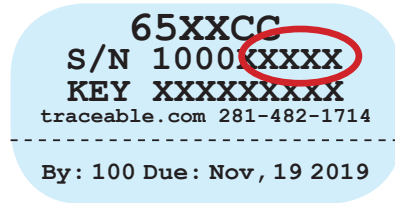

## Open Web Browser on your **MOBILE DEVICE**

8

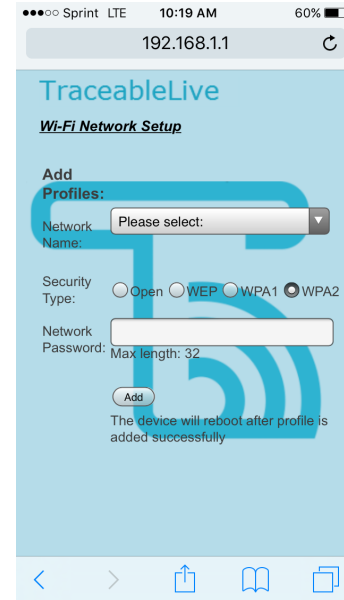

Enter into web browser: **192.168.1.1**


Select desired local network to which the device  
will be connected, and enter password for network.

The most common network is **WPA2**.

## Press Add/Confirm device restart

9

Click Add

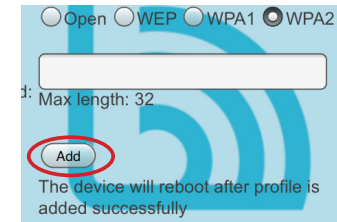

Unit will reboot, connect to WiFi, and begin posting data.

**Note:** Upon successful connection, you will receive an error  
message in your browser (Example: Safari cannot open  
the page because the network connection was lost.), and  
device reboot may take up to 1 minute, mobile will exit WiFi.

**Note:** The WiFi symbol will flash on device display until  
the device posts data successfully.

**Note:** ALLOW UP TO 30 MINS FOR DEVICE DATA TO  
UPDATE AND START TRANSMITTING DATA.

If and ERR message occurs on device, press acknowl-  
edge button (✓) and repeat steps 4-7.

Log-in to TraceableLIVE account to manage device  
settings, alarms, users, and payment options.

Device and account setup is now complete.

92-6500-30

TraceableLIVE®  
QUICK START  
GUIDE

# Create Account

1

Visit <https://www.traceablelive.com>  
Click on New Account.

# Enter Account/User Information

2

Fill in all required fields.  
**Note:** Password must be a minimum of 8 characters including upper and lower case letters, numbers and special character (!@#%\$%^&\*).

Click Register

Register

# Enter Payment Info./ Confirm Charge

3

If this screen does not appear, go to log in screen, account settings and add payment information.  
**Note:** Payment must be set up prior to adding a device to your service.


Click Save Payment Info

# Enter Device Serial Number Info (located on back of unit)

4

Go to [www.traceablelive.com](http://www.traceablelive.com)  
Click on Add New Device

Enter S/N & Device Key  
Click SUBMIT

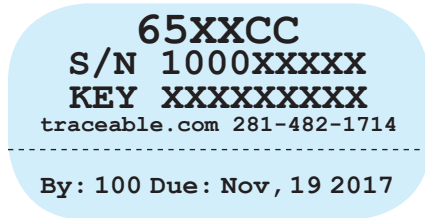

S/N = Serial Number  
KEY = Device Key

# Enter Location, and Alarm Information

5

Complete all required fields.  
Click Add